

Things to do before you leave:

- On the Linux workstation close the FE Beam Permit
- Click the ABSB Open Allowed to No
- Call Control Room x8899 and ask them to close the Port Shutter
- Save all your files (copy them to the network drive, P, and to your USB stick), exit the OLIS program and shut down the computer
- Switch off the OLIS System CD controller box (next to the PC) and all other equipments (temperature controller Quantum box, Peltier pump etc).
- Close the N₂ gas valve to a standby at 5L/min.
- Switch off computer monitors (Linux WS, and the ring status display as well) and all other electronic equipments (Nanodrop, PCs in DP room, etc.)
- Switch off lights by the main door.
- Check that everything is done.